



Updating your AAA software to the latest revision

DO NOT uninstall the old version of AAA


1. Double click on the AAA setup installer for the latest revision. DO NOT run as administrator.

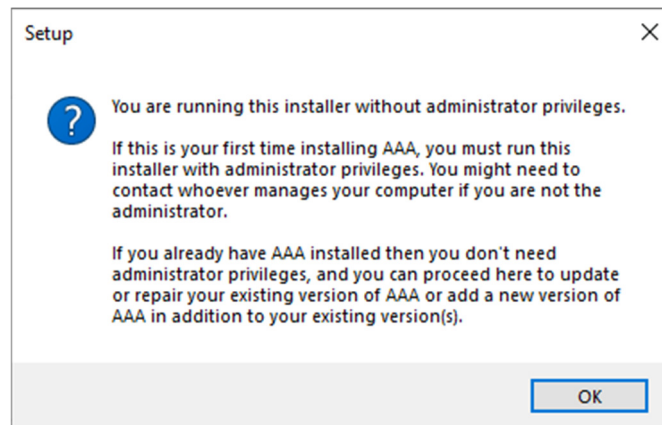


2. Select I accept the agreement and Click **Next**

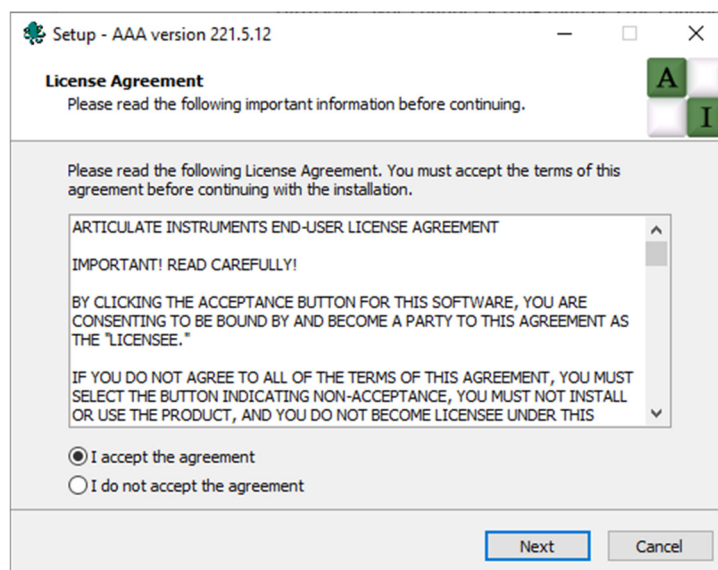


3. Click **Next** unless you have reason to install the new revision elsewhere. This will overwrite the old version.

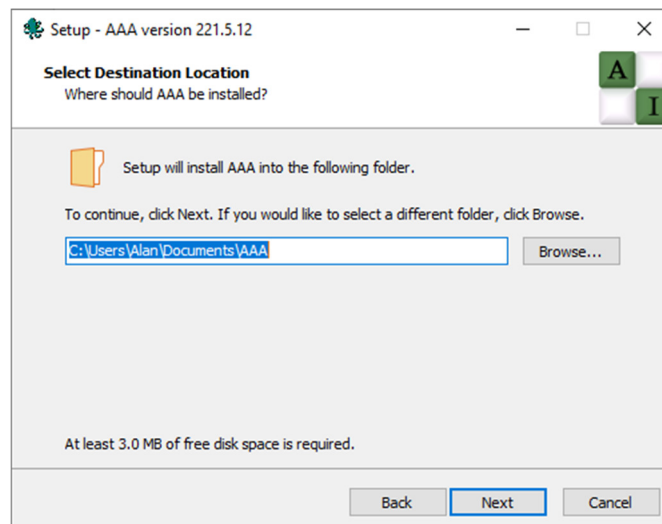


4. Click Yes unless you have reason to install the new revision elsewhere. This will overwrite the old version.

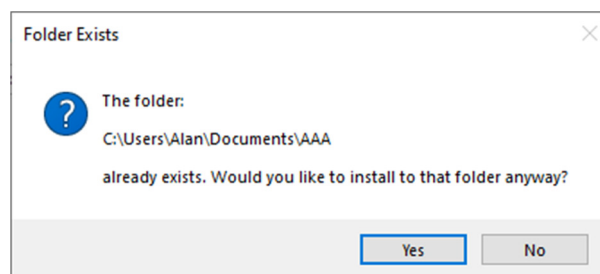


5. Click Next to install the revised version of AAA, up-to-date tongue and lip models, and default AAA settings for this revised version.

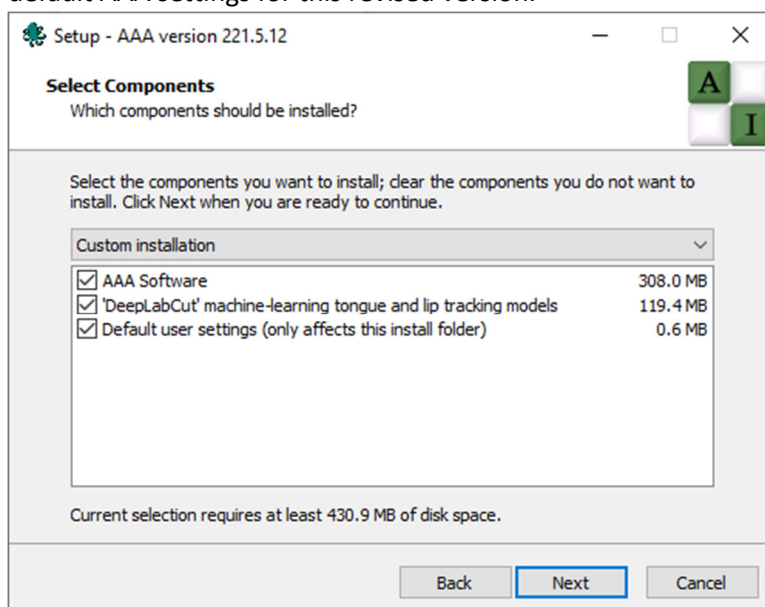

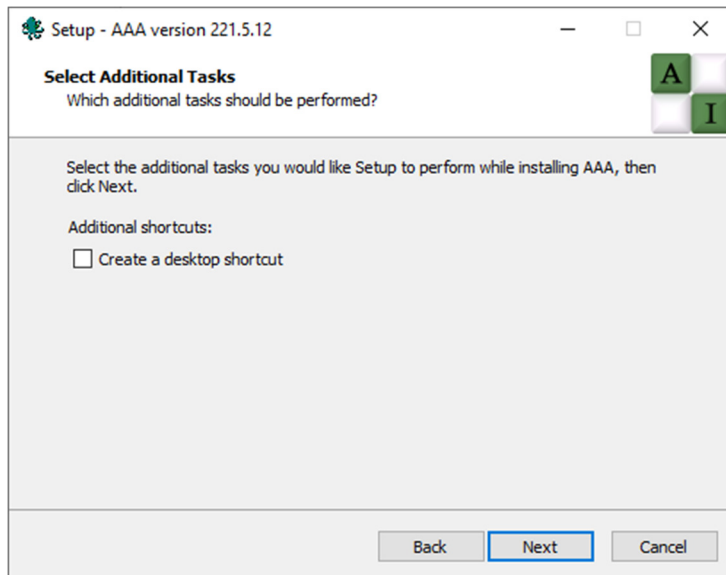


6.

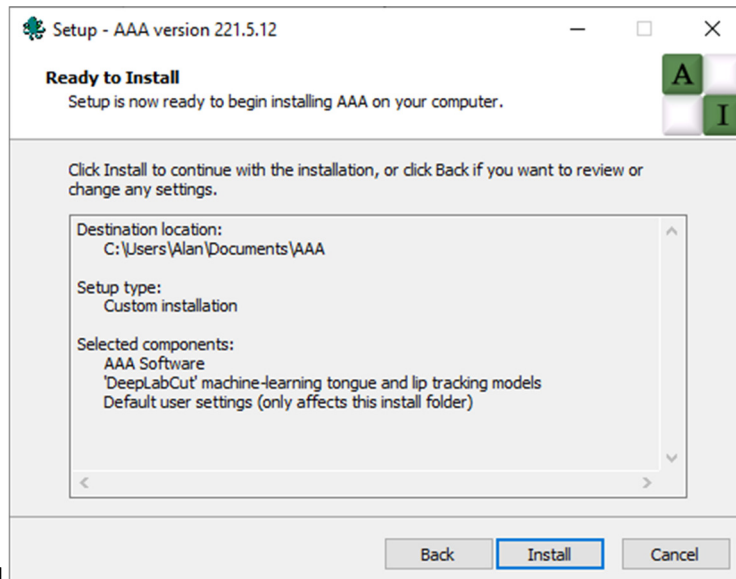


7. Click Install

8. CLICK CANCEL as Miniconda has already been installed.

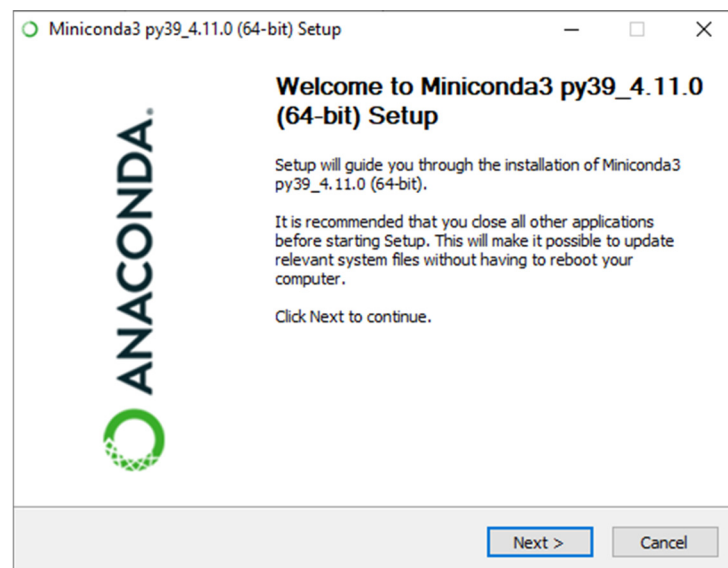

9. AAA will automatically start to check that DeepLabCut is installed.

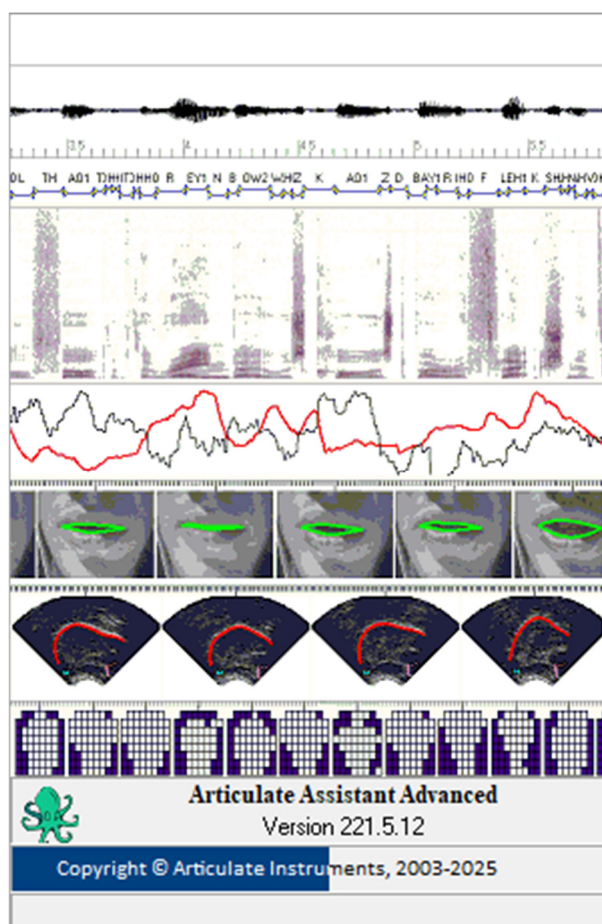

Click

