

Focusrite Scarlett soundcard configuration

Unfortunately, Focusrite change their Scarlett model designs every couple of years, so there are slightly different instructions for connecting audio depending on which model you have.

The following settings apply to all models:

1. The 48V phantom power button must be switched ON or there will be no microphone signal.
2. The channel that the PStretch is connected to must be set to LINE and not INST or AIR
3. The monitor dial must be set to the 2 o'clock setting to provide the right speaker level.



4. The Direct monitor must be OFF.
5. Speaker or headphones must be connected to the LINE OUTPUTS on the rear of the soundcard.
6. Scarlett must be selected in Windows sound manager for both playback as well as recording.



Focusrite Scarlett Solo Gen2



Focusrite Scarlett Solo Gen3



Focusrite Scarlett Solo Gen4



Focusrite Scarlett 4i4 Gen3



Focusrite Scarlett 4i4 Gen4

Differences:

1. For all except the Focusrite Scarlett Solo Gen2 two 6.25mm-phono/RCA adapters should be plugged into the LINE OUTPUTS. This allows the twin phono/RCA to 3.5mm jack cable to be connected.



1. For all except Scarlett Solo Gen4, the microphone should be connected to channel 1.